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Version 1.3 09/2023

User instructions Microscope camera

Before use

You should ensure that the device is not exposed to direct sunlight, temperatures which are too high or too low, vibrations, dust or a high level of humidity. The ideal temperature range is between 0 and 40°C and a relative humidity of 85% should not be exceeded.

Always make sure that you use an approved power cable. Thus, possible damages by reason of the development of overheating (fire hazard) or an electric shock can be prevented.

Do not open the housing and touch the internal component. There is the risk of damaging them and affecting the functionality of the camera.

In order to carry out cleanings always disconnect the power cable from the camera.

Always keep the sensor clear from dust and do not touch it. Otherwise, there is the risk of affecting the microscopic image. In case of non-use always attach the protective covers.

Technical data

Model KERN	Resolution	Interface	Sensor	Frame rate	Colour / Mono- chrome	Supported operating systems
ODC 825	5,1 MP	USB 2.0	1/2,5" CMOS	6,8 -55 fps	Colour	Win, XP, Vista, 7, 8, 10
ODC 831	3,1 MP	USB 3.0	1/3" CMOS	27,3 – 53,3 fps	Colour	Win, XP, Vista, 7, 8, 10
ODC 832	5,1 MP	USB 3.0	1/2,5" CMOS	14,2 – 101,2 fps	Colour	Win, XP, Vista, 7, 8, 10
ODC 841	20 MP	USB 3.0	1" CMOS	15 - 60 fps	Colour	Win, XP, Vista, 7, 8, 10

Scope of delivery

- Microscope camera
- USB cable
- Object micrometre for calibration
- Software CD

Free download:

www.kern-sohn.com > DOWNLOADS > SOFTWARE > Microscope VIS Pro

- Eyepiece adapter (Ø 23,2 mm)
- Adjustment rings (Ø 30,0 mm + Ø 30,5 mm) for eyepiece adapter

Mounting

1. Remove the black cover at the bottom of the camera.
2. The thread, where the cover was attached, is a standardised C mount thread. Thus, there are special C mount adapters needed for the connection to a microscope.
3. For the mounting to the microscope the C mount adapter is attached to the connection point of the microscope. After that the camera has to be screwed onto the C mount adapter.

Important:

The choosing of the right C mount adapter depends on the used microscope model. It must be an adapter, which is adjusted to the construction of the microscope and recommended by the manufacturer as appropriate for the relevant microscope.

4. If necessary, adjust the microscope according to the trinocular usage (with help of the trino toggle rod / trino toggle wheel).

PC connection

1. Establish USB connection via USB cable.
2. Installing the software with help of the CD / Download.
3. The software-internal "User Guide" includes all information and instructions about the operation of the software or of digital microscopy.

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